

CRIAC QDI 2010

Microspectrophotometer

1) Operator Safety:

- a. Take care when loading/unloading samples to insure hands and fingers are clear prior to moving the stage.
- b. Do not touch lamp housing, as it gets hot.
- c. Wear UV glasses when using it.
- d. Report safety concerns to the NFCF staff through e-mail or by calling 3-8001.

2) Avoiding Machine Damage:

- a. Follow procedures carefully.
- b. Careful not to damage optics when using microscope.
- c. Use care when loading and unloading samples from holders to avoid damage.
- d. Report any system contamination to the NFCF staff.

3) Potential System Hazards:

- a. Pinch Points
- b. Compressed gasses.
- c. Ultraviolet Light.
- d. AC Voltage.
- e. Hot Surfaces.

For Emergency Assistance call 412-624-2121. (Internal 4-2121)